

In vitro Toxicology

Structural 3D cardiotoxicity assay

Background Information



Numerous studies have shown that cell responses to drugs in 3D culture are improved from those in 2D, with respect to modeling *in vivo* tissue functionality, which highlights the advantages of using 3D-based models for preclinical drug screens'

⁶Nam KH, Smith AS, Lone S, Kwon S and Kim DH (2015) *J Lab Autom*; In press

- Drug induced cardiovascular toxicity is the leading cause of attrition during drug development. Drugs can exert functional toxicities such as arrhythmia and morphological (structural) damage to the myocardium¹. Evaluation of the potential for both types of cardiotoxicity by novel compounds is essential for the discovery of safe drugs.
- The myocardial tissue comprises 30% cardiomyocytes and 70% non-myocytes, the majority of which are endothelial and fibroblast cells. These non-myocytes are essential to myocardial structure and function^{2,3} with emerging evidence suggesting important roles within drug induced cardiovascular toxicity⁴.
- Mitochondrial disruption, calcium dyshomeostasis and cellular ATP content have been identified as major targets for structural cardiotoxins⁵.
- Three dimensional (3D) confocal HCS allows the simultaneous detection of each cell health parameter in combination with a measure of cellular ATP.

Protocol

Microtissue

Human induced pluripotent stem cell derived cardiomyocytes (iPSC-CM's), cardiac endothelial cells and cardiac fibroblasts

Analysis Platform

Confocal Cellomics ArrayScan® XTI (Thermo Scientific)

Test Article Concentrations

8 point dose response curve with top concentration based on $100x C_{max}$ or solubility limit. 3 replicates per concentration.

Test Article Requirements

50 μL of a DMSO solution at a concentration of 200x top concentration (top concentration = 100x C_{max}) or equivalent amount in solid compound

Time Points

72 hours (Others available on request)

Quality Controls

Negative control: 0.5% DMSO (vehicle) Positive controls: Sunitinib (Ca²⁺ homeostasis) and dasatinib (mitochondrial membrane potential)

Data Delivery

Minimum effective concentration (MEC) and AC_{50} value for each measured parameter (microtissue count, microtissue size, DNA structure, calcium homeostasis (Ca^{2+}), mitochondrial mass (Mito Mass), mitochondrial membrane potential (MMP) and cellular ATP content)

Figure 1

Representative 3D confocal high content screening (HCS) images of isoproterenol calcium dyshomeostasis in spontaneously beating cardiac 3D microtissues labelled with Hoechst (blue) to detect DNA structure, Fluo-4 AM (green) to detect calcium dyshomeostasis and TMRE (red) to detect mitochondrial disruption.











Compound	C _{max} (μΜ)	<i>In vivo</i> toxicity	hESC-CM prediction (Pointon et al 2013) ⁵	H9c2 monolayer	H9c2 MTs	iPSC-CM's monolayer	Tri-cultured cardiac MTs	Most sensitive
				MEC (µM)				feature
Dasatinib ⁷	0.72	Structural cardiotoxin Non- structural cardiotoxin	Positive structural cardiotoxin	0.529	NR	7.32	2.08	MMP
Doxorubicin ⁸	15.34			0.04	0.115	0.04	0.04	ATP
Fluorouracil ⁹	4.61			1.88	NR	1.4	0.0407	Ca ²⁺
Idarubicin HCI10	0.12			<0.04	<0.04	0.04	<0.04	ATP
Imatinib Mesylate11	3.54			13.7	3.53	29.3	22.6	ATP
Lapatinib ⁷	4.18			4.57	8.33	0.04	5.9	ATP
Sunitinib Malate12	0.25			0.896	0.114	0.04	0.817	Ca ²⁺
Cyclophosphamide13	153.20		Negative structural cardiotoxin	NR	NR	NR	30.8	Mito Mass
Isoproterenol HCl ¹⁴	0.01			NR	NR	NR	2.1	Ca ²⁺
Acyclovir	6.66			NR	NR	NR	NR	-
Buspirone HCI	0.03			NR	0.237	NR	NR	Microtissue size

Table 1

Structural cardiovascular toxicity prediction of 12 reference compounds categorised according to literature data.

Tri-culture cardiac 3D microtissues (MTs), H9c2 3D microtissues (MTs) and H9c2 monolayers were incubated with test compound for 72 hours. The cell models were analysed using the confocal mode of Cellomics ArrayScan® XTI (Thermo Scientific) following which cellular ATP content was measured using CellTiterGlo® (Promega).

 $\label{eq:MEC} \begin{array}{l} \mathsf{MEC} = \mathsf{minimum} \ \mathsf{effective} \ \mathsf{concentration} \\ \mathsf{NR} = \mathsf{no} \ \mathsf{response} \end{array}$

Figure 2

Graphical representation of (a) sunitinib induced calcium dyshomeostasis and (b) dasatinib induced mitochondrial membrane potential disruption in spontaneously beating cardiac 3D microtissues.

All reference compound toxicities were correctly predicted in the spontaneously beating cardiac tri-culture 3D microtissue model including isoproterenol (MEC 2.1µM, calcium dyshomeostasis (Table 1 and Figure 2)) and cyclophosphamide (MEC 30.8µM, mitochondrial mass (Table 1)) which previously went undetected by Pointon *et al* (2013)⁵ and Cyprotex's in-house H9c2 data.

Control compound sunitinib displays cytosolic calcium increase (calcium dyshomeostasis) followed by gross cytotoxicity (microtissue loss) (Figure 2a) while control compound dasatinib displays mitochondrial membrane potential loss without gross cytotoxicity (microtissue loss) (Figure 2b).



The combination of an *in vitro* 3D model that better recapitulates the *in vivo* cellular physiology of the myocardium with a multiparametric HCS and cytotoxicity assay presents a viable screening strategy for the accurate detection of novel therapeutics that cause drug induced structural cardiovascular toxicity early in drug development.

References

- ¹ Laverty HG, et al., (2011) Br J Pharmacol **163(4)**; 675-693
- ² Brutsaert DL, (2003) Phys Revs 83; 59-115
- ³ Souders CA, et al., (2009) Circ Res **105**; 1164-1176
- ⁴ Mikaelian I, et al., (2010) Tox Sci **117(1)**; 144-151
- ⁵ Pointon A, et al., (2013) Tox Sci **132(2)**; 317-326
 ⁶ Nam KH, et al., (2015) J Lab Autom; In press
- ⁷ Force T, et al., (2007) Nat Rev Cancer 7; 332-344

- ⁸ Minotti G, et al., (2004) Methods Enzymol **378**; 340-361
- ⁹ Schimmel KJ, et al., (2004) Cancer Treat Rev 30(2); 181-191
- ¹⁰ Anderlini P, et al., (1995) J Clin Oncol **13(11)**; 2827-2834
- ¹¹ Kerkelä R, *et al.*, (2006) *Nat Med* **12(8)**; 908-916
- ¹² Chu TF, et al., (2007) Lancet **370**; 2011-2019
 ¹³ Floyd JD, et al., (2005) J Clin Oncol **23(30)**; 7685-7696
- ¹⁴ Zhang J, *et al.*, (2008) *Toxicol Pathol* **36(2)**; 277-278